

Maximo Dry PCR Master Mix**Cat. S295; 12x8-Tube Strips***Ready-to-use lyophilized PCR-Master-Mix in 8-tubes-strips***Features:**

- All components like Taq polymerase, dNTP's, reaction buffers, enhancers and stabilizers are lyophilized
- Transportation and at room-temperature
- simple, and fast setting up procedure for high yield and repeatable results

Description:

Maximo Dry Master-Mix is optimized and ready-to-use mixture of all components for a successful PCR. Only your primers and your DNA Template has to be added to the 20 µl assay.

Maximo Dry Master-Mix contains a thermostable DNA polymerase that possesses a 5'→3' polymerase activity and a double-stranded specific 5'→3' exonuclease activity. The enzyme consists of a single polypeptide with a molecular weight of 94kDa.

List of components:

12x8 lyophilized PCR-Tubes (0,2ml) "flat cap",

Unit definition:

One unit incorporates 10 nmol of deoxyribonucleotide into acid-precipitation material in 30min at 74°C

Quality control:

- PCR with various templates – genomic DNA, Phage Lambda DNA
- 2 kb DNA amplification from 50 ng DNA
- batch variation and level of bacterial DNA contamination

Transportation: at room temperature

Storage: store at room-temperature in an aluminium coated bag or on another dry place; humidity < 65 % when sealing is opened

Usage:

Components	Volume per reaction
1 Tube of Maximo Dry Mastermix	
Forward Primer:	0,2-1 µM – 0,4-2µl / 10 µM
Reverse Primer	0,2-1 µM – 0,4-2µl / 10 µM
Template DNA	1-50 ng
Sterile dest. Water (molecular grade)	up to 20 µl total reaction volume

Note:

- vortex all solutions and spin down carefully before using

GeneON .. a good decision ..

Contact Phone +49-(0)-621- 5720 864 Fax: +49-(0)-621-5724 462

E-Mail: <mailto:info@geneon.net> WEB: <http://www.GeneOn.net> Version: 10.2009 UK

Unless specified otherwise, all products of GeneON are sold for research use only.

.. a good decision ..

General Thermo-Cycler protocol:

Step	Time	Temperature
Initial denaturation	2 min	94°C
30 Cycles:		
Denaturation	25-30 sec	94°C
Annealing ^{x1}	25-30 sec	50-68°C
Elongation ^{x2}	0,5-3 min	72°C per 1kb
Final extension	2 min	70-72°C

Note:

- ^{x1} depends on the melting temperature of the primers
- ^{x2} depends on the length of the fragments (1 min/kb)

Ordering information

Cat.-no	Description	Amount
S295	Maximo Dry Master Mix	12x8 Flat Cap PCR-tubes

.. a good decision ..

GeneON .. a good decision ..

Contact Phone +49-(0)-621- 5720 864 Fax: +49-(0)-621-5724 462

E-Mail: <mailto:info@geneon.net> WEB: <http://www.GeneOn.net> Version: 10.2009 UK

Unless specified otherwise, all products of GeneON are sold for research use only.